



## Quick Guide: Nirvascan ASP-NIR-M-Reflect ( internal #NIR-S-G1)

### Product Overview



#### STEP 1 : Turn on the NIR-S-G1

Move the switch to the right side, the green and blue LEDs will light up indicating that the device and Bluetooth are on. The scanner is ready to use.



#### STEP 2 : Start your scanning experience

Turn on your mobile device App and connect NIR-S-G1 via Bluetooth; or PC (with SW) connect NIR-S-G1 via USB. Put the object close to the scan window, then press the button. Within a few seconds, you will be able to view the scan results on your App or SW.



#### Remark A : Charge NIR-S-G1

When the charge LED is blinking, please plug the Micro-B USB cable to get the device charged from any standard USB port or USB charger.



#### Remark B : Disable Bluetooth function

If you don't need to use the Bluetooth function, you can disable it by pressing the button for at least 3 seconds, when you release the button the blue LED light will turn off.

